

The final date for return of results is indicated on the request/report form dispatched with the samples.

Contact details:

The Organisers - FEPTU
HPA Microbiology Services Colindale
61 Colindale Avenue,
London, NW9 5EQ, UK.
Fax: +44 (0) 20 8200 8264
Tel: +44 (0) 20 8327 7119
e-mail: foodeqa@hpa.org.uk
www.hpa.org.uk/eqa/standard



Results should be returned on-line using the web-based reporting system. Go to: www.kpmd.co.uk/hpa/. You will need your log-in details for this process. Please contact us if you experience problems with the on-line system.

Standard Scheme - Instruction Sheet

Each distribution consists of **two** evacuated glass vials containing mixtures of freeze-dried micro-organisms. The vial contents require reconstitution prior to examination.

Refer also to the Safety Data Sheet: www.hpa.org.uk/eqa/standard

Storage:

Store the samples at **2 - 8°C** on receipt.

Safety Recommendation:

The samples contain micro-organisms of ACDP Hazard Groups 1 and 2. Your laboratory must undertake a risk assessment before handling these samples with particular consideration to the conditions and procedures for opening evacuated vials to prevent the risk of aerosol.

Opening Vials:

Remove the plastic tear-off seal by positioning the vial with the arrow on the plastic flip-top pointing away from you, flip the top up then carefully but deliberately pull downwards and to the right (or the left if you are left handed) until the seal separates; then still holding onto the plastic top, gently remove altogether and discard into a sharps container.

Reconstitution:

The samples are reconstituted in **200 ml** nutrient broth (**pre-warmed** to 30 ± 2 °C) as follows:

- Aseptically remove the rubber stopper from the vial of the freeze-dried sample.
- Transfer approximately 1 ml of the **200 ml** warmed nutrient broth to the vial and allow to stand for 2 - 5 minutes.
- Transfer the vial contents back to the remaining nutrient broth.
- Rinse the vial a further 3 - 4 times with the **200 ml** nutrient broth sample.
- Shake the reconstituted samples before examination.

Examination:

- **Each reconstituted sample is equivalent to a 200 g food sample.**
- Undertake the examinations listed on the request/report form within 45 minutes of reconstitution.
- Examine in accordance with routine procedures.

Reporting:

- Return your results on-line using the web-based reporting system. Go to: www.kpmd.co.uk/hpa/
- You will need your log-in details for this process.
- Contact us if you experience problems with the on-line system.

Return of Results:

- The distribution will be closed on the deadline date for return of results.
- Results cannot be accepted after the distribution has closed.